



Bee Stinger

March 2012

INSTRUCTION MANUAL

Gastight Syringe for the BASi Bee Syringe Pump

2701 Kent Avenue
West Lafayette
Indiana 47906

www.BASiInc.com

Bee Stinger

A Gastight Syringe for the BASi Bee Syringe Pump
Model-00001

Warning

Laboratory syringe for research purposes. Not for human use.

Important! Before using this syringe remove the plunger from the barrel. Dip plunger into deionized or distilled water and re-insert into glass barrel. Move the plunger up and down five times. Movement of the syringe plunger in a dry barrel may cause excessive wear on the Teflon seat, reducing the lifetime of the seal and causing leaks.

Cleaning: Deionized or distilled water, acetone or methyl alcohol are all recommended cleaning solvents for this syringe. Avoid detergents, phosphates and alkalines whenever possible. Syringes may be soaked overnight in a mild solvent or immersed in an ultrasonic cleaner for a few hours if necessary